

# GPC Filters

Positively Charged, Single Layer  
PES Membrane




Fine Particle Removal

Endotoxin Removal

*GPC filters are part of Critical Process Filtration's economical, general service product line that can be used to lower the total cost of filtration from simple to the most complex applications. Produced using the same quality materials and manufacturing excellence as our other product lines, you can be assured of their performance, dependability and scalability. While general service filters are not integrity tested or validated for retention, these highly efficient filters remove large amounts of contaminants early in the process to reduce the load on your expensive downstream filters. They are also useful as stand-alone clarifying or particle removal filters in less critical applications.*

GPC cartridge and capsule filters are made with a single layer positively charged Polyethersulfone (PES) membrane. They are used for the filtration of aqueous fluids that contain negatively charged biological contaminants, such as endotoxin, virus and other cell fragments. GPC filters are well suited for endotoxin and fine particle retention, delivering high flow and throughput. Pore sizes range from 0.03 to 1.2  $\mu\text{m}$ .

Critical Process provides unrivaled delivery times, technical consulting before purchasing, and very competitively priced high-performance products. Our comprehensive testing & analysis and validation services support your team whenever they need it. Your process experts partnering with our filtration experts is how we deliver your company's solution right the first time.

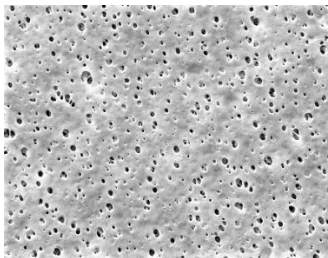

GPC filters are recommended for endotoxin and fine particle retention in:

- Process Water
- DI Water



CARTRIDGES – Nominal Dimensions

Length: 5 to 40 in. (12.7 to 101.6 cm)

Outside Diameter: 2.75 in. (7.0 cm)



CAPSULES – Nominal Dimensions

Length: 2 to 30 in. (5.1 to 76.2 cm)

Outside Diameter: 3.50 in. (8.9 cm)

## Maximum Operating Parameters

	CARTRIDGES	CAPSULES
<b>Liquid Operational Pressure</b>	N/A	80 psi at 68 °F (5.52 bard at 20 °C)
<b>Gases Operational Pressure</b>	N/A	60 psi at 68 °F (4.14 bar at 20 °C)
<b>Operating Temperature (water)</b>	180 °F at 30 psid (82 °C at 2.07 bard)	110 °F at 30 psid (43 °C at 2.07 bard)
<b>Forward Differential Pressure</b>	80 psid at 68 °F (5.52 bard at 20 °C) (Liquid and Gas)	Liquid - 80 psid at 68 °F (5.52 bard at 20 °C) Gas - 60 psi at 68 °F (4.14 bar at 20 °C)
<b>Reverse Differential Pressure</b>	50 psid at 68 °F (3.45 bard at 20 °C)	50 psid at 68 °F (3.45 bard at 20 °C)
<b>Recommended Changeout Pressure</b>	35 psid (2.41 bard)	35 psid (2.41 bard)

## Sanitization & Sterilization

<b>Filtered Hot Water*</b>	90 °C (194 °F), 30 minutes, multiple cycles, max 3 psid forward flow	N/A
<b>Inline Steam*</b>	275 °F (135 °C), 30 min, 25+ cycles	N/A
<b>Autoclave*</b>	250 °F (121 °C), 30 min, 25+ cycles	250 °F (121 °C), 30 min, 25+ cycles
<b>Chemical Sanitization</b>	Performed using industry standard concentrations of hydrogen peroxide, peracetic acid, sodium hypochlorite and other selected chemicals.	

\*Cartridge Filters – For all elevated temperature procedures above, a stainless-steel support ring is required.

## Filtration Area (Nominal)

	CAPSULES	CARTRIDGES AND CAPSULES				CARTRIDGES
		2"	5"	10"	20"	
Length	5.08cm	12.7cm	25.4cm	50.8cm	76.2cm	101.6cm
Area	1.2 ft <sup>2</sup>	3.4 ft <sup>2</sup>	7.3 ft <sup>2</sup>	14.6 ft <sup>2</sup>	21.9 ft <sup>2</sup>	29.2 ft <sup>2</sup>
	0.11m <sup>2</sup>	0.32m <sup>2</sup>	0.68m <sup>2</sup>	1.36m <sup>2</sup>	2.04m <sup>2</sup>	2.72m <sup>2</sup>

## Construction Materials

<b>Filtration Media</b>	Single Layered Asymmetric Positively Charged Polyethersulfone (PES) Membrane
<b>Media Support</b>	Polypropylene
<b>End Caps, Center Core, Outer Support Cage, Capsule Housing</b>	Polypropylene
<b>Sealing Method</b>	Thermal Bonding
<b>O-Rings/Gaskets Cartridges only</b>	Buna, Viton® (or FKM), EPDM, Silicone, FEP Encapsulated Silicone, FEP Encapsulated Viton (or FKM)

## Extractables

GPC filters typically exhibit low levels of non-volatile residues.



## Non-Fiber Releasing

The GPC filters comply with Title 21 CFR sections 210.3 (b)(6) and 211.72, for non-fiber releasing filters.

## FDA and EC Compliance



Materials meet the requirements listed by the FDA as appropriate for use in articles intended for repeated food contact as specified in Title 21 CFR sections 174.5, 177.1500, 177.1520, 177.1630, 177.2440, and 177.2600 as applicable. All materials used to make the filters are listed in European Commission Regulation EU/10/2011, Annex 1.

## Flow Rates for GPC Cartridges by Pore Size

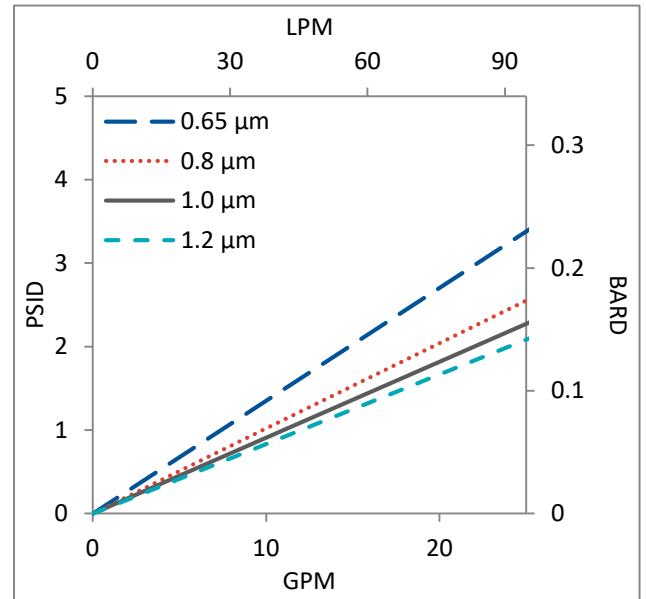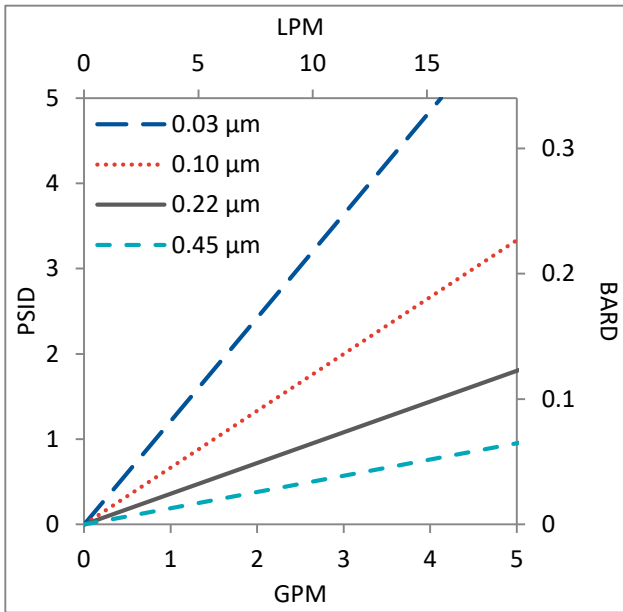

Flow rates for Cartridge filters are per 10-inch length. The test fluid is water at ambient temperature.

## Flow Rates for GPC Capsules by Pore Size

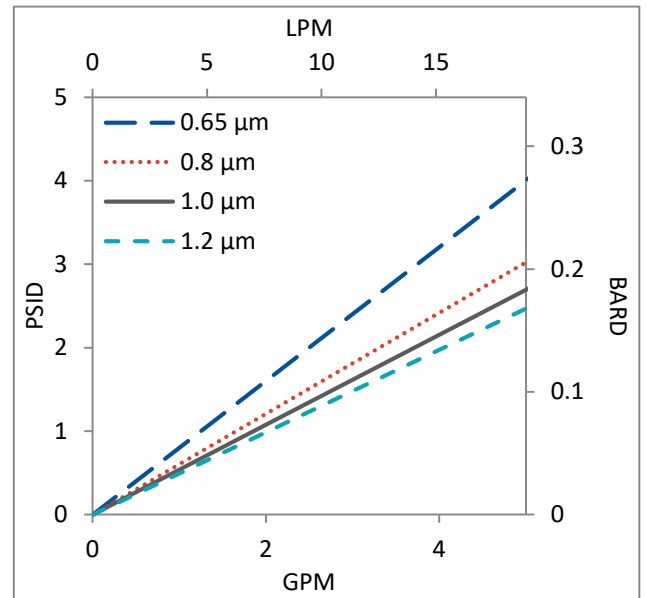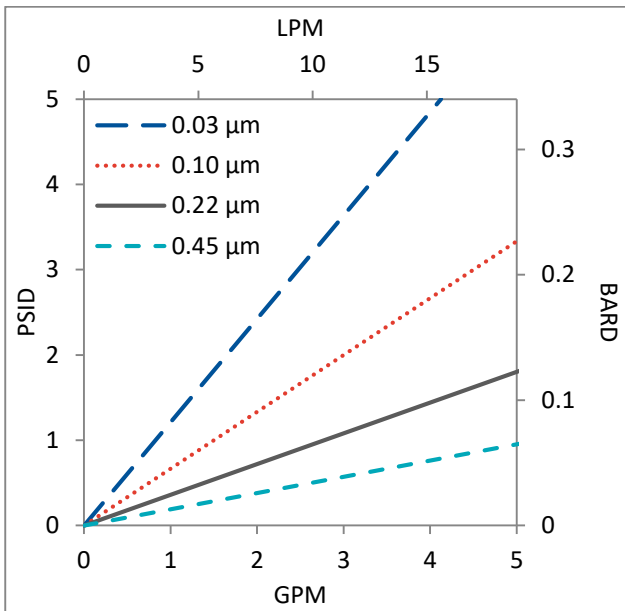


Flow rates for Capsule filters are tested using a 2" capsule filter with 1" sanitary inlet and outlet ports. The test fluid is water at ambient temperature. Flow rates for larger capsules will scale with filtration area. Rates will vary based on end configuration of the capsule.

## GPC Filters Ordering Information


All Critical Process filters are configurable to meet customer specifications.  
Fill in the corresponding codes in the boxes below to build your Part Number.

To consult with one of our technical team members, request a quote or place an order:  
call (603) 880-4420 Ext. 106, or send an email to [sales@criticalprocess.com](mailto:sales@criticalprocess.com)

### Cartridge Filters

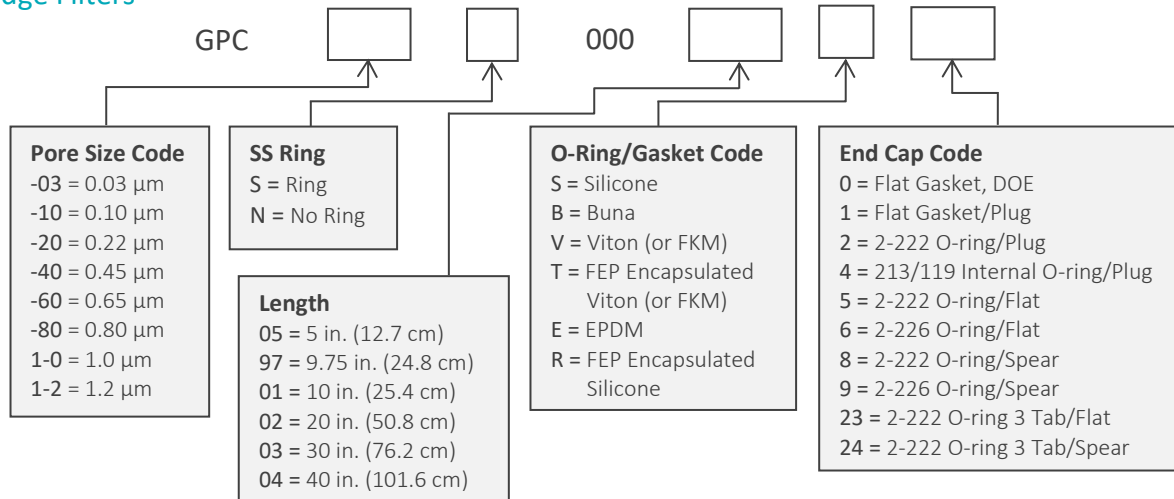

### Capsule Filters

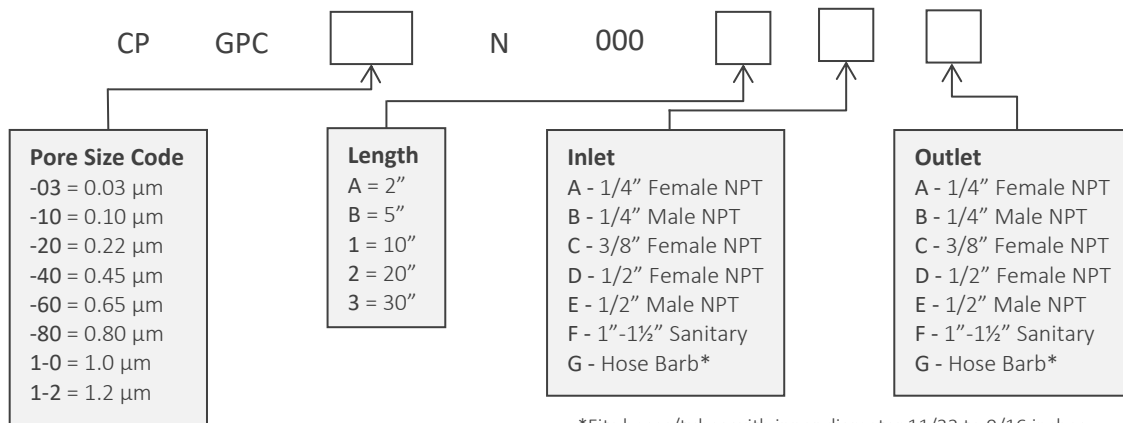

### Housings

CPF offers Model CSH sanitary housings and Model CIH industrial housings in Single-Round (Inline and T-Style) and Multi-Round configurations. Contact us or visit our website for more information.



One Chestnut Street  
Nashua, NH 03060  
603.880.4420  
FAX: 603.880.4536  
CriticalProcess.com

The information contained herein is subject to change without notice. The Critical Process Filtration logo is a trademark of Critical Process Filtration, Inc. Viton is a trademark of DuPont Performance Elastomers L.L.C.  
© 2021 Critical Process Filtration, Inc. • All Rights Reserved

Data Sheet GPCDS Rev B